




ALARM BUZZER

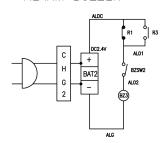



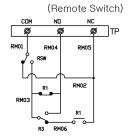

lf the Auto Back-up System is Installed (LN2)

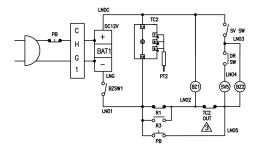

BZ2	DC12V Buzzer				
R1,3	Support Relay				
TC20UT	Temperature Controller Output				
DRSW	DRSW DOOR Switch				
PT2	PT2 RTD Sensor (Pt100Ω)				
TC2	TC2 Temperature Controller				
SV5	Solenoid Valve (LN2)				
SVSW	SVSW Solenoid Valve Switch				
PB					
BZ1	DC12V Buzzer				
BZSW1	Buzzer Switch(AL-1)				
BAT1	Battery(DC12V)				
CHG1	Battery Charger				
SIMBOL					
If the Auto Back-up System is Installed		(LN2)			
R1,3	Support Relay				
BZ3	DC 3V Buzzer				
BZSW2	BZSW2 Buzzer Switch(AL-1)				
BAT2					
CHG2	Battery Charger				
SIMBOL DESCRIPTION		REMARKS			
Alarm					
R1.3					
RSW					
TP	TP Terminal,Output Signal				
SIMBOL	SIMBOL DESCRIPTION				
	Remote				
THSW1-2	THSW1-2 Thermo Start Switch				
TR					
SV1-4					
PSW1-2	PSW1-2 Pressure Control Switch				
RL Error Lamp(AL2)					
T2,3 Timer					
T1	T1 Compressor Delay Timer				
COMP1-2					
TH1-2					
F1SB.F2SB	SB.F2SB Support Fan Motor				
F1-2	F1-2 Fan Motor				
CS					
R1-5					
MG1-2	MG1-2 Magnet Relay				
PT1	, ,				
TC1					
PSW	PSW Power Switch				
SIMBOL	SIMBOL DESCRIPTION				

Description: Wiring Diagram CRYO 170/230



Drawing no.:	5086084	Revision:	05
Drawing (initials):	DBL	Date:	201210
Approval (initials):	MET	Date:	201210

All rights to this drawing belong to Arctiko A/S. The drawing may not be copied nor presented to any other person without prior written consent of Arctiko A/S.