

ACCPTFRZ



PerformaTherm™ Medical Transport Cooler Replacement Freezer PCM Panels

Six (6) -21°C Freezer PCM Panels

This manual outlines the procedure for use of the ACCPTFRZ -21°C. It includes the conditioning of the -21°C PCM Panels for -25°C to -15°C transport and provides step by step assembly instructions. As it is a universal pack-out, conditional and assembly are the same for both Summer and Winter conditions and are as follows:

CONDITIONING


1. Ensure all PCM-21°C panels are frozen by storing them in a freezer at $-30^{\circ}\text{C} \pm 3^{\circ}\text{C}$ for 24 Hours or until the panels are clearly frozen solid.
2. Immediately proceed to pack out procedure

PACK OUT PROCEDURE

1. Place one (1) PCM-21°C panel at the base of the VIP tub.
2. Place one (1) PCM-21°C panel at each of the sides of the VIP tub.
3. Place products inside the payload created in steps 1 & 2.
4. If needed, add dunnage to fill the void space.
5. Place one (1) PCM-21°C panel on top of the four PCM-21°C panels. It will fit into the beveled edges of the other PCM-21°C panels to sit flush.
6. Fully zip the top of the PerformaTherm™ ACCPT3L or ACCPT5L to ensure it is completely closed.

Optional: Use a digital data logger to monitor the payload for accurate temperature sensitive product(s) and regulatory compliance.

PACK OUT DIAGRAM

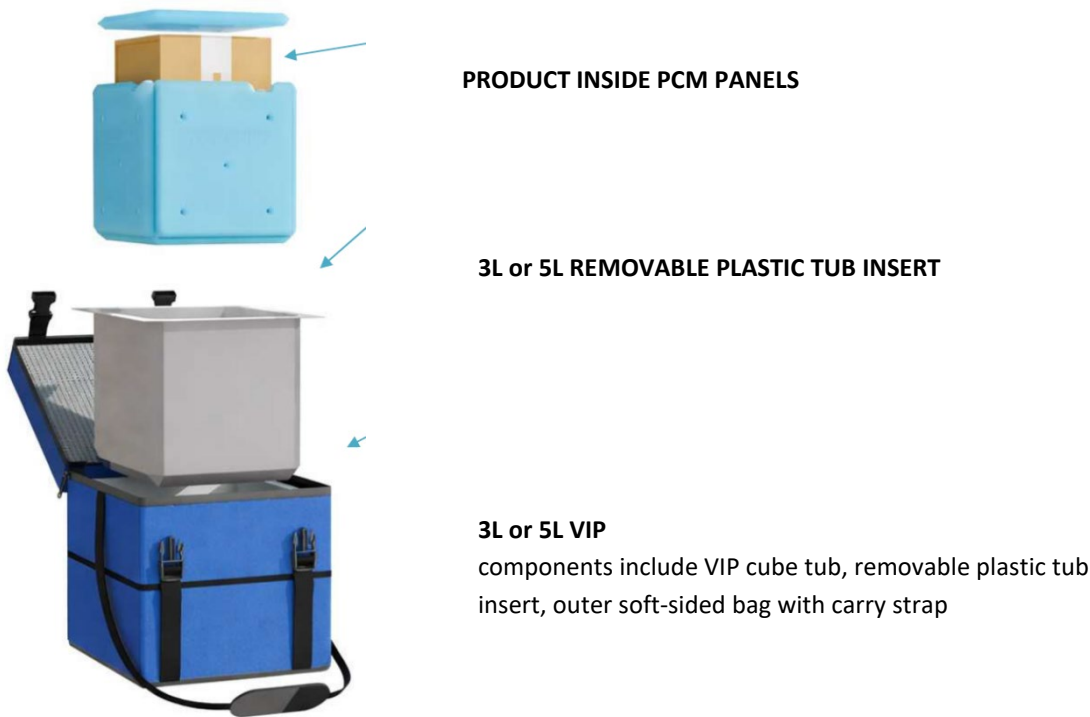

Exterior Cooler Dimensions	3L capacity: 11.5" x 11.5" x 10.5" 5L capacity: 14.5" x 11.5" x 10.5"
Payload Dimensions	3L 5"x5"x5" 5L 9"x5"x5"
Validated Performance	ISTA 7E Summer/Winter performance tested with Vacuum Insulated Panel (VIP) Temp-Shield™ system. Maintains accurate and stable temperature (-21°C) for 2+ days.
Applications	Blood, plasma & organ transport Vaccines & biologics Clinical trial materials Hospital, laboratory & field operations
Limited Warranty	PERFORMA-THERM™, 1-year Limited Warranty – covering defects in materials and workmanship from delivery date. Exclusion: not valid if altered, misused, damaged by accident, or due to negligence. To make a warranty claim, the purchaser must contact Thermco at info@thermcoproducts.com , or by calling 1-877-900-4141.
Extended Warranty	Additional one (1) year from date of purchase. Contact Thermco Products for additional details.

HTS Code 3926.90.9989 Manufactured Country of Origin – China

SDS PCM+5 panel.pdf