



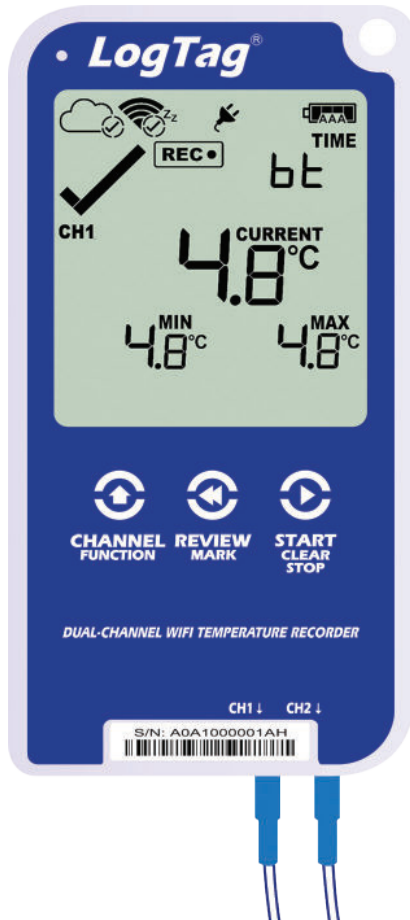

UTRED30-WIFI

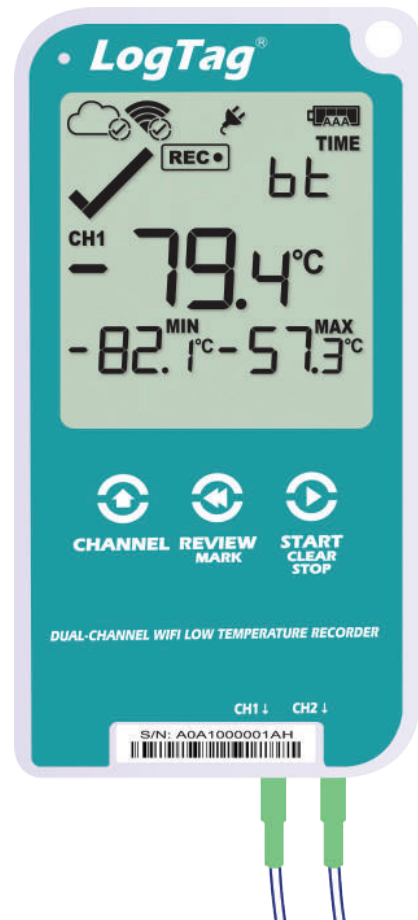

UTREL30-WIFI

LogTag®

Bluetooth® Configuration Guide

Version B - August 2022

www.logtag.com

Registering the Device

The LogTag Mobile application allows users access to LogTag Online using a mobile device. Through Bluetooth® capabilities the application allows your device to be connected to the network and be remotely configured. Profile and timezone can be configured remotely for the logger.

For more information or alternative configuration methods refer to your Wi-Fi loggers quick start guide or user manual on their product pages.

Downloading the LogTag Mobile Application

1. Navigate to either the App Store or Google Play store within your mobile device.
2. Search for the LogTag Mobile application.
3. Download the latest LogTag Mobile application.




4. Launch the application and follow the sign in process.

Ensure your device is either plugged in to your computer or an AC adapter using the USB cable provided.

Note: Allow all permissions for the applications use of location, Bluetooth® and camera.

Registering the Device (Continued)

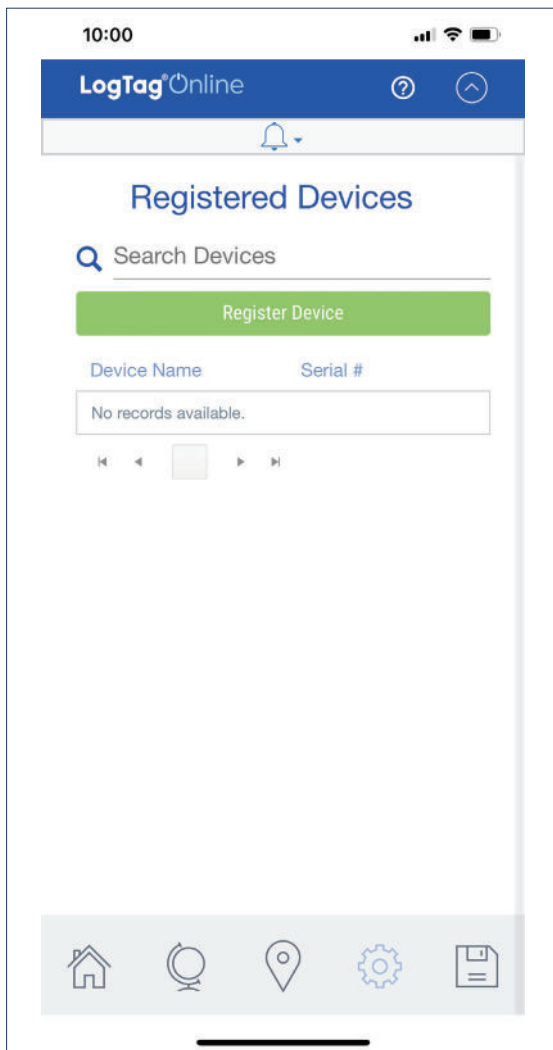


Ensure you have a location slot available on your LogTag Online account.

5. Navigate to the devices tab  and select 'Register Device'.

6. Scan or enter the barcode of your LogTag logger, then select Next.

Note: Hold the CHANNEL FUNCTION button to activate Bluetooth® if 'bt' does not show.

Registering the Device (Continued)

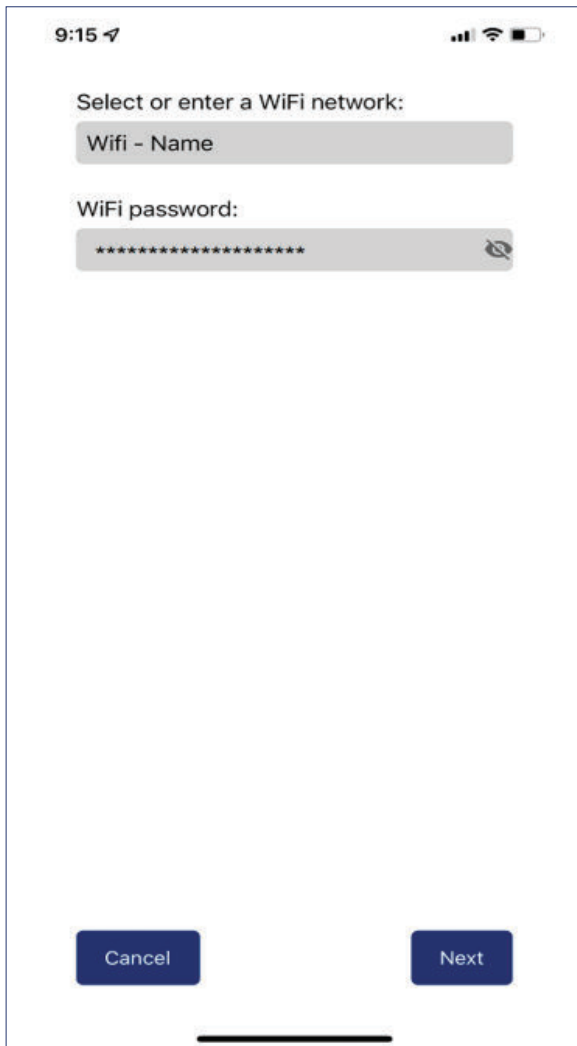


9:15

Select or enter a WiFi network:

Wifi - Name

WiFi password:

Cancel Next

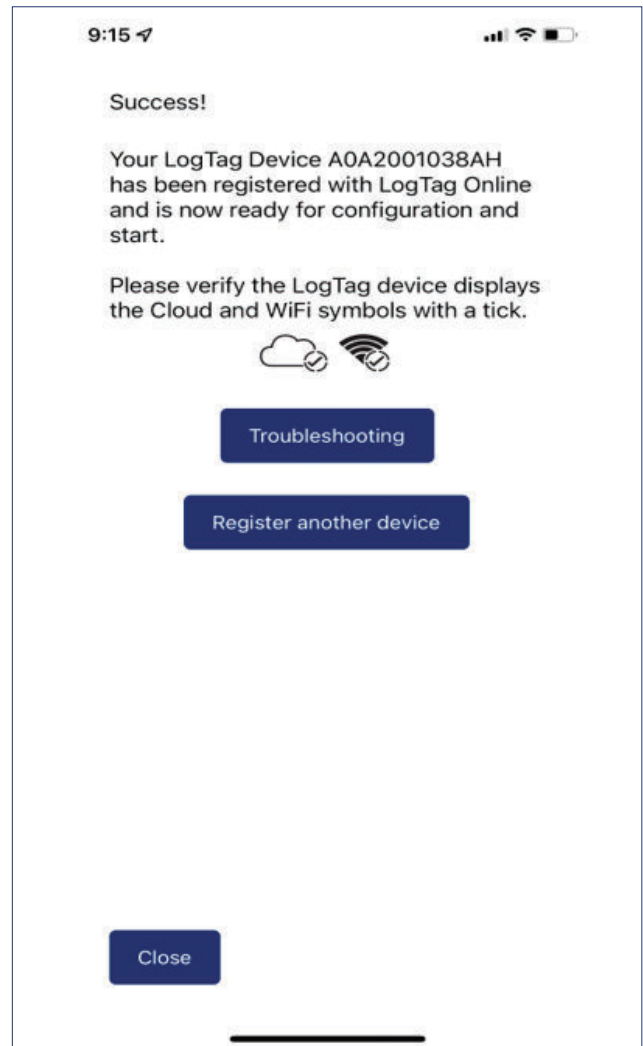

9:15

Success!

Your LogTag Device A0A2001038AH has been registered with LogTag Online and is now ready for configuration and start.

Please verify the LogTag device displays the Cloud and WiFi symbols with a tick.

Troubleshooting

Register another device

Close

7. Connect your device to the network by entering or selecting your Wi-Fi network name and providing the network password, then select 'Next'.

Note: Ensure the Wi-Fi password is correct. To view the password select the eye icon.

8. Your device will connect to your local network using your provided Wi-Fi credentials, once completed select 'Close'.

Select 'Troubleshooting' if you are experiencing issues.